

Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922) 49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Ижевск (3412)26-03-58
Иваново (4932)77-34-06
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Киргизия (996)312-96-26-47

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Россия (495)268-04-70

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сыктывкар (8212)25-95-17
Сургут (3462)77-98-35
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Казахстан (772)734-952-31

Тольяти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Улан-Удэ (3012)59-97-51
Ульяновск (8422)24-23-59
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

<https://grant.nt-rt.ru> || gtq@nt-rt.ru

УФ-камера для ПЦР UVT-B-AR



The Grant UVT-B-AR is an economical general purpose PCR UV cabinet in for carrying out short-term procedures.

The UVT-B-AR cabinets are equipped with an open UV lamp installed in the upper part of the hood. UV-radiation from these lamps effectively disinfects the working area chamber, inactivating DNA / RNA fragments during 15-30 minutes of exposure. The UV light switches off automatically when door is opened. The model has a stainless steel work surface, metal framework and shock proof glass panels.

The built-in UV cleaner-recirculator both increases the maximum density of UV light and generates 25 m³/h air flow exchange – prevents unwanted contamination and protects the user from direct UV light during manipulation. The UV cleaner-recirculator consists of UV lamp, dust filters and a fan. The recirculation effect increases the maximum density of the UV light making it sufficiently effective for DNA / RNA inactivation. The recirculator processes 360 UV-cabinet volumes per hour, creating permanently aseptic conditions within the cabinet.

The digital timer controls the duration of the direct UV irradiation. An additional daylight lamp provides a bright, clear working area.

- 1 x 25W 254 nm open UV lamp
- High intensity UV cleaner – 25m³/hour cleaner-recirculator continuous air flow with 1cm UV irradiation distance
- UV exposure controlled by 24h digital timer
- Convenient, easy to use digital timer for accurate control of UV exposure
- White lamp provides local illumination of the workplace to optimise visual control during operations
- Automatic switch-off of irradiated UV light when door opened
- Supplied with 1 access port and internal power socket

Applications

Life science – germicidal and virucidal, inhibition of DNA and RNA contamination, applications requiring no residual decontaminants such as disinfectants, operations with DNA/RNA amplicons, microbial research.

Specifications

	UVT-B-AR	UVT-B-AR INL
Weight	32 kg	32 kg
Construction	stainless steel frame and working area	stainless steel frame and working area
Panels	glass with UV-protective film	glass with UV-protective film
Front opening with three adjustable positions	•	•
Open UV lamp, 25 W bactericidal, 254 nm, ozone free	1	1
Open UV lamp, 30 W bactericidal, ozone free	–	–
Bactericidal air recirculator, 25 m ³ /h air flow exchange	•	•
UV recirculator, 25W (efficiency >99% per 1 cycle)	1	1
UV recirculator, 30W (efficiency >99% per 1 cycle)	–	–

White lamp for workplace illumination 15W	1	1
White lamp for workplace illumination 30W	—	—
Digital timer (non-stop) 0 to 24 hours	•	•
Internal power outlets	1	—
Internal working area	645mm x 490mm	645mm x 490mm
Flow rate m3/h	7	7
Nominal operating voltage V	120 or 230 (50-60Hz)	120 or 230 (50-60Hz)
Power consumption W	67 (0.3A)	67 (0.3A)

Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922) 49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Ижевск (3412)26-03-58
Иваново (4932)77-34-06
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Киргизия (996)312-96-26-47

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Россия (495)268-04-70

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сыктывкар (8212)25-95-17
Сургут (3462)77-98-35
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Казахстан (772)734-952-31

Тольяти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Улан-Удэ (3012)59-97-51
Ульяновск (8422)24-23-59
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

<https://grant.nt-rt.ru> || gtq@nt-rt.ru